**Stereomikroskop
Stereo Microscope
Stéréomicroscope
Estereomicroscopio**

**Bedienungsanleitung
User Manual
Mode d'emploi
Instrucciones de uso**

**Modell / Model
Modèle / Modelo 33208**

ESCHENBACH

Deutsch	3
English.....	11
Français	19
Español.....	27

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Index

Safety instructions	11
Proper use	13
Items supplied	13
Warranty	13
The Operating Elements of the Microscope	14
Working with LED Ring illumination.....	15
Transmitted light operation.....	15
Taking the microscope into use	16
Maintenance and care	16
Disposal	17

Safety instructions

Risk of electric shock

- Connect the illumination units of the appliance exclusively via its power adapter to a correctly installed mains power socket supplying a voltage of 230 V~ / 50 Hz.
- Do not operate the illumination units if the power adapter, the adapter's power cable or the appliance itself is damaged. Check the power adapter from time to time for signs of damage. Arrange for a defective power adapter to be repaired or replaced as soon as possible by Customer Services.
- You must not open or repair the adapter housing yourself. There are no user-serviceable elements inside. Should you do so, appliance safety can no longer be assured and the warranty will become void.

- Protect the appliance from moisture and humidity. Use it only in dry rooms, do not use it outdoors or close to liquids.
- Make certain that neither the appliance nor the mains adapter could become wet or damp or sustain any other damage during use.
- Notice regarding separation from mains-power
To completely separate the appliance from mains power, the power adapter must be unplugged from the wall socket. For this reason, the appliance should be placed where unhindered access to the power socket is assured at all times, so that in an emergency you can immediately withdraw the plug. To eliminate the risk of fire, prevent unintended activation and save energy costs, the power adapter should be unplugged from the wall socket when the appliance is not in use.
- Always grasp the mains adapter directly when you wish to unplug it. Do not pull on the cable and never touch the mains adapter with wet hands, this could result in a short circuit or you receiving an electric shock. Ensure that the cable cannot be trapped in any way. Never tie knots in the power cable, and do not join it to other cables. Place the power cable where it cannot be trodden on.

 ***Risk of accidents!***

- Do not allow children to handle optical appliances without supervision, as they cannot always assess the potential injury hazards.
- Ensure that no one can stumble over the illumination cable.
- Should the illumination become damaged, the appliance may no longer be taken into use. Arrange for the appliance to be checked and, if necessary, repaired by qualified technicians.
- This appliance is not intended for use by individuals (including children) with restricted physical, physiological or intellectual abilities or deficiencies in experience and/or knowledge unless they are supervised by a person responsible for their safety or receive from this person instruction in how the appliance is to be used. Children should be supervised to ensure that they do not play with the appliance.

Risk of suffocation!

- Do not leave packaging material lying unattended. Plastic foils and bags and pieces of styropor etc. are not toys, they can be dangerous for small children.

Proper use

This appliance is intended exclusively for use as a microscope. No warranty claims will be granted for damage resulting from improper use!

Do not subject the appliance to high temperatures, direct sunlight, severe vibrations, shocks or tremors or high mechanical loads.

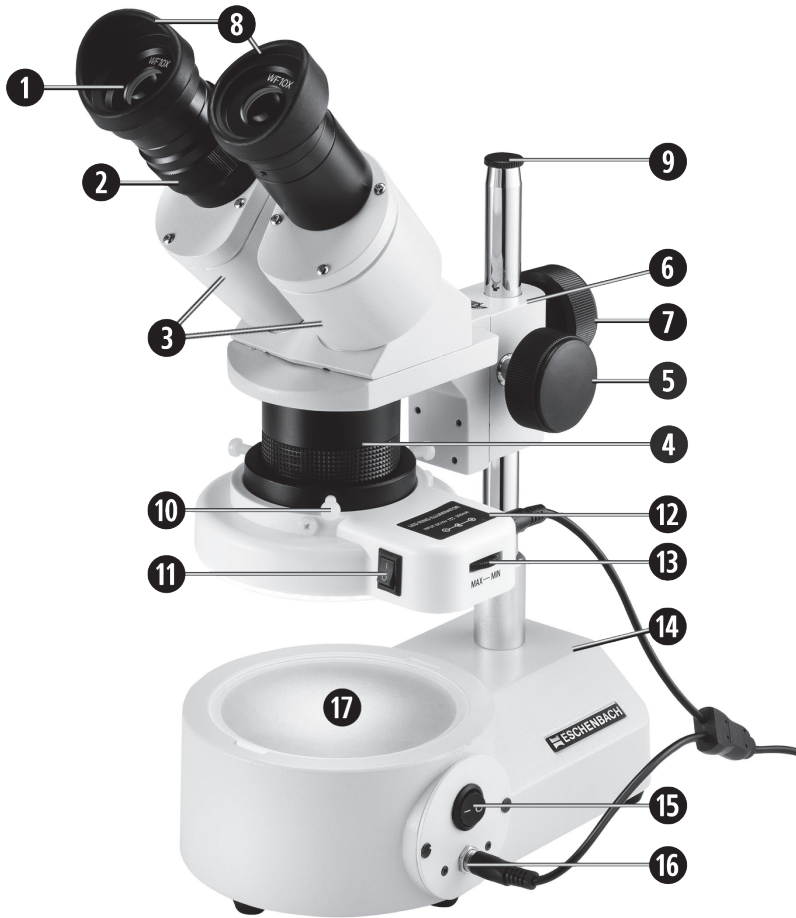
Items supplied

- ▶ Microscope head complete with focussing gear box and transmitted light illumination
- ▶ LED Ring illumination for reflected light operation
- ▶ Mains power adapter
- ▶ Table stand
- ▶ Object bearer disc, made of glass, for transmitted light operation
- ▶ Dust protection cover
- ▶ 2 Eyepiece cups
- ▶ Hook wrench
- ▶ Operating instructions

Warranty

We vouchsafe within the scope of current statutory regulations the function of the appliance described in these instructions in regard to defects which may occur due to manufacturing or material flaws. We accept no liability for damage occasioned by improper handling, falls or impacts of all kinds. Warranty claims must be supported by the sales receipt!

The Operating Elements of the Microscope


- ① Eyepieces
- ② Dioptic adjuster
- ③ Eyepiece bearer
- ④ Lens barrel
- ⑤ Gear wheel for height adjustment
- ⑥ Gear box with stand receiver
- ⑦ Securing wheel (for locking the microscope head)
- ⑧ Eyepiece cups
- ⑨ Cap screw
- ⑩ Plastic screw
- ⑪ On/off switch for LED ring illumination
- ⑫ LED ring illumination
- ⑬ Dimmer for LED ring illumination
- ⑭ Table stand
- ⑮ On/off switch for transmitted light illumination
- ⑯ Power supply socket
- ⑰ Object bearer disc

Working with LED Ring illumination

Amongst the items supplied is an LED Ring illumination unit 12 for illuminating the objects being examined by microscope. To use this, proceed as follows:

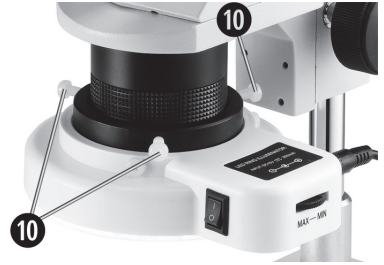- Unscrew the 3 plastic screws 10 out of the LED Ring illumination 12 until they no longer project in the opening for the lens barrel 4.
- As shown on the right, push the ring illumination over the lens barrel and then secure it by carefully tightening the 3 plastic screws.
- Insert the power adapter into a correctly installed mains power socket and then connect one of the plugs of the adapter cable with the socket on the ring illumination.
- Place the On/Off switch 11 of the LED Ring illumination at the position '1' to switch the appliance on in reflected light operation.
- Adjust the required light output with the MAX-MIN regulator 15.
- After use, switch the LED Ring illumination off by placing the On/Off switch at the '0' position.

Transmitted light operation

- For transmitted light operation use the supplied object bearer disc 17 made of glass.
- Insert the power adapter into a properly installed mains power socket. Insert one of the plugs of the adapter cable into the power supply socket 16.
- Place the On/Off switch 15 of the illumination unit at the position '1' to switch the appliance on in transmitted light operation.
- After use, switch the illumination unit off by placing the On/Off switch at the position '0'.

Both illumination units use the same power supply via a Y cable.
Both illumination units can be used simultaneously.

Taking the microscope into use

The appliance is delivered largely pre-assembled.

- Place the two eyepiece cups ③ onto the eyepieces ①.
- Unscrew the protective cap of the lens barrel ④ anti-clockwise.
- Look through the two eyepieces. Adjust the eyepiece spacing to your eye spacing by turning the eyepiece bearer ③ either inwards or outwards.
- For the individual optimal contrast select either the black or the white side of the object bearer disc ⑰.
- Place the object that is to be examined by microscope on the object bearer disc.
- With largish objects that are to be observed or processed, the spacing-adjustment using the lateral gear wheels ⑤ may not be sufficient. In this case, you should first of all carry out a rough distance adjustment by using the securing wheel ⑦. For this, orientate yourself on the following table:

Objective magnification	Eyepiece magnification	Total magnification	Work spacing [mm]	Field of view Ø [mm]	Flat image field = Field of view Ø [mm] x Objective magnification
2 x	W 5 x	10 x	75	10,5	21
4 x	W 5 x	20 x	45	5,5	22
2 x	W 10 x	20 x	75	10	20
4 x	W 10 x	40 x	45	5	20
2 x	W 15 x	30 x	64	7	14
4 x	W 15 x	60 x	43	3,5	14
2 x	W 20 x	40 x	64	4,5	9
4 x	W 20 x	80 x	43	2,5	10

W = Broad field

The 10 x Eyepiece is supplied, the others are optionally available.

- Select the required magnification (2x or 4x) by turning the lens barrel through 180° for each.
- For fine adjustment of the focus, use the gear wheel ⑤ to adjust the distance of the microscope head to the object, this will ensure an optimally sharp image.
- Modify the dioptic adjuster ② to meet your personal requirement.

Maintenance and care

- For cleaning, do NOT use aggressive chemical substances, solvents or abrasives, these could damage the surfaces.
- To avoid soiling of the appliance, make use of the supplied dust protection cover when it is not in use.
- Should the appliance nonetheless become soiled, clean it with a soft and lightly moistened cloth.
- For cleaning the eyepieces, use a lens cleaning cloth.

- Exercise care when cleaning the eyepieces, the upper surfaces can be easily scratched.

By using the supplied hook wrench you can adjust the free movement of the gearing. On one side of the gear wheel axis there is a small bushing with corresponding holes. Make the required adjustment by turning it either clockwise or anti-clockwise.

Disposal

Under no circumstances should the appliance be disposed of with domestic waste at the end of its useful life. This product is subject to the provisions of European Directive 2012/19/EU.

Arrange for the product, or parts of it, to be disposed of by a professional waste disposal company or by your communal waste facility. Observe the currently applicable regulations. In case of doubt, please contact your waste disposal centre. Dispose of packaging materials in an environmentally responsible manner.

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Eschenbach Optik GmbH